



# Bacterial and Fungal Whole Genome Re-sequencing

## Service Introduction

Bacterial and fungal whole genome re-sequencing is a critical tool to complete the genomes of known bacterium and fungi, as well as to compare multiple genomes or to map genomes of new organisms. It is of great importance to sequence entire genomes of bacterium and fungi in order to generate accurate reference genomes, to do microbial identification and other comparative genome studies.

Compared with traditional methods like PCR, next generation sequencing (NGS) excels in the way of multiplexing vast number of samples, providing identifications of SNP/InDel/SV with great reliability, turnaround time and cost performance.

## Service Specifications



### Library preparation and sequencing strategy

- Genomic DNA or PCR product will be fragmented to prepare 350 bp insert library
- Paired-end (PE) 150 bp strategy by using Illumina NovaSeq platform



### Quality standard and sequencing depth


- No less than 80% of bases with a Q30 (99.9% accuracy) quality score
- Recommend sequencing coverage of  $\geq 100\times$  for bacterial genomes,  $\geq 50\times$  for fungal genomes



### Turnaround time and analysis options

- Typically 18 business days from confirmation of library preparation to raw data delivery
- Various analysis options from standard bioinformatics package to custom analysis

## Project Workflow

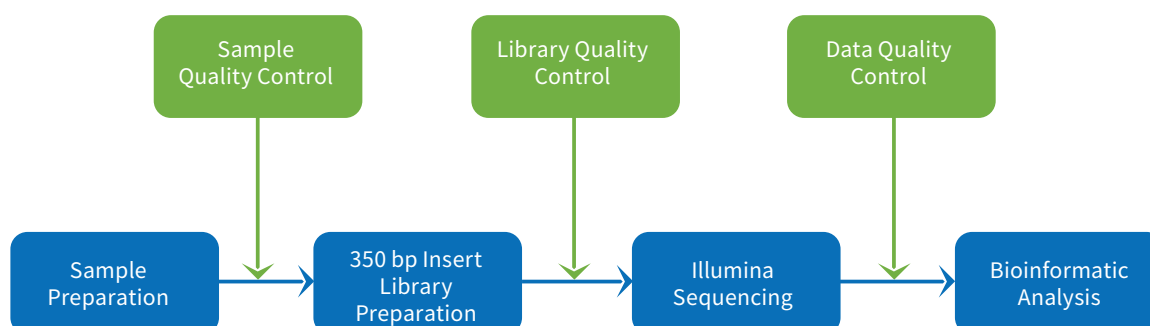


## Sample Requirements

Sample Type	Amount	Volume	Concentration	Purity
Genomic DNA	≥ 300 ng	≥ 20 µL	≥ 5 ng/µL	OD260/280=1.8-2.0, No degradation or RNA contamination
PCR Product	≥ 1.5 µg	≥ 20 µL	≥ 20 ng/µL	OD260/280=1.8-2.0, No degradation or RNA contamination

## Data Analysis

Besides raw data output, Novogene offers a variety of standard and customized analysis for your bacterial and fungal genome re-sequencing projects.

### Standard Analysis Pipeline

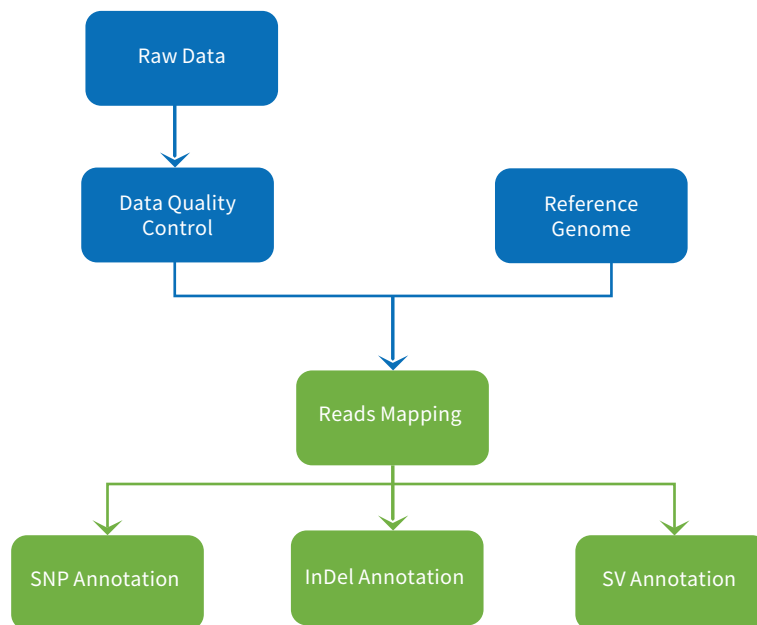

- Data quality control: e.g. filtering reads containing adapter or with low quality
- Alignment with reference genome, statistics of sequencing depth and coverage
- SNP/InDel/SV calling, annotation and statistics

### Custom Analysis Options

A wide range of customized bioinformatics analysis are available, please contact your Novogene representative.



Follow us on LinkedIn

## Novogene Corporation Inc.

📍 8801 Folsom Blvd #290, Sacramento, CA 95826

☎ 916-252-0068-383   ✉ [inquiry\\_us@novogene.com](mailto:inquiry_us@novogene.com)   🌐 [en.novogene.com](http://en.novogene.com)

For Research Use Only. Exclusive for Clients in North and South America.

Copyright©2011-2021 Novogene Corporation. All Rights Reserved.

Information and specifications are subject to change at any time without notice:

Please contact your Novogene representative.